

# Bio-Tape™

## Standardized Method for Surface Sampling

### Sampling Procedure


Use this line as a guide to let you know where the middle of the adhesive area is.



Remove PEEL liner from adhesive area and discard. Do not replace after sampling.

Document the sample location here.

1. Remove slide from provided slide mailer
2. Document the sample location on the slide label and complete any required documentation for the laboratory.
3. Peel off protective liner from slide to expose adhesive
4. Place slide with the adhesive side down on the surface to be sampled. (Use the centering line to orient the adhesive over the exact sample location desired)
5. Gently press down to ensure contact is made. DO NOT press down hard. The adhesive is very tacky and only needs to be pressed gently.
6. Carefully remove slide from surface and place back into slide mailer. DO NOT replace the plastic "Peel" liner.
7. Send slide to laboratory for analysis.

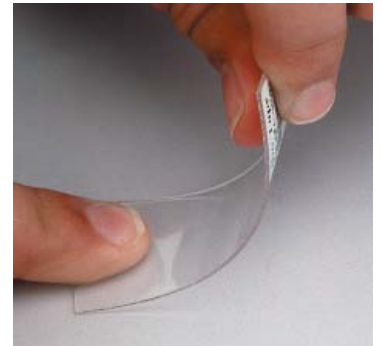

Flexible plastic slide allows you to place the adhesive on the surface to be sampled while holding the non-sticky end of the slide.

#### Note:

It is recommended that multiple samples are collected from both suspect and non-suspect (control) areas for comparison.



Sampling Equipment Specialists

1-800-282-0073  
2860 23rd Avenue North  
St. Petersburg, FL 33713  
www.zefon.com